



SECTION A-A
SCALE 1 : 4

The Duraline Recirculating Lab Faucet comes with the Recirculating Module and Tube temporarily assembled.

- STEP 1.** Remove (unscrew) the Recirculating Module and Recirculating Tube from the through deck nipple.
- STEP 2.** Mount the faucet in the desired location through a 1" diameter hole. Tighten to deck with the Ring Nut supplied.
- STEP 3.** Insert the Recirculating Tube into and through the Through Deck Nipple until the 3/8" male NPT thread mates with the 3/8" female NPT thread in the inlet connection of the Through Deck Nipple. Tighten by hand plus 1/4 wrench turn.
- STEP 4.** Connect a 3/8" supply line tube to the 3/8" inlet compression tube fitting.
- STEP 5.** Connect a 1/4" return line tube to the 1/4" outlet compression tube fitting.
- STEP 6.** Pressure test all connections.

Note: A high density PTFE tape is the recommended thread sealant.

UNLESS OTHERWISE SPECIFIED DIMENSIONS ARE IN INCHES		2950 Airway Avenue, Costa Mesa, CA 92626		SCALE: N/A	Format: SD	SHEET 1 OF 1
---	--	--	--	------------	------------	--------------

Designer		TITLE: Duraline Recirculating Lab Faucet, Deck Mount, With Recirculating Module
Review		
Engineer		

	DWG. NO.	DL-DDR-2 (Natural Polypropylene) DL-DDR-3 (Natural PVDF)	REV
--	----------	---	-----

5 4 3 2 1